

TS 097 – AMPICILLIN SUPPLEMENT

INTENDED USE

For isolation of *Aeromonas hydrophila* from foods.

COMPOSITION

Ingredients	Concentration
Ampicillin	10mg

(Per vial sufficient for 500 ml medium)

APPLIED IN MEDIUM

TM 1074 - SA AGAR BASE (as per APHA)


TM 1137 - AEROMONAS STARCH DNA AGAR BASE

INSTRUCTION FOR USE

Rehydrate the contents of 1 vial aseptically with 2 ml of 50% v/v alcohol. Mix well and aseptically add to 500 ml of sterile, molten, cooled (45-50°C) TM 1136 - Aeromonas Isolation Medium Base. Mix well and pour into sterile petri plates.

STORAGE:

Vials should be stored in sealed container at 2-8°C.

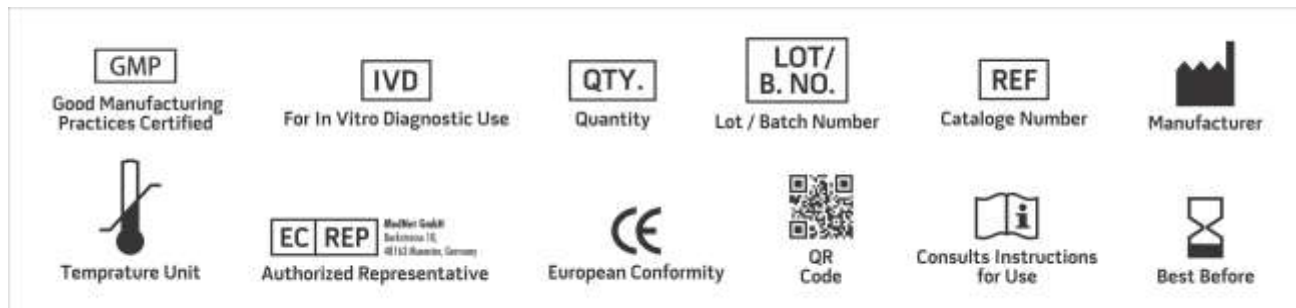

NOTE: Please consult the Material Safety Data Sheet for information regarding hazards and safe handling Practices.

***For Lab Use Only**
Revision: 02 Feb., 2022