

TS 111 – GEORGE KIMMIG SELECTIVE SUPPLEMENT

INTENDED USE

For selective isolation and cultivation of fungi.

COMPOSITION

Ingredients	Concentration
Penicillin G	20000IU
Streptomycin	20mg

(per vial sufficient for 500 ml medium)

Applied in Medium


TM 1211 - KIMMIG FUNGI AGAR BASE

INSTRUCTION FOR USE

Rehydrate the contents of 1 vial aseptically with 5 ml of sterile distilled water. Mix well and aseptically add it to 500 ml of sterile, molten, cooled (45-50°C) TM 1211 - KIMMIG FUNGI AGAR BASE. Mix well and pour into sterile petri plates.

STORAGE

Vials should be stored in sealed container at 2-8°C.

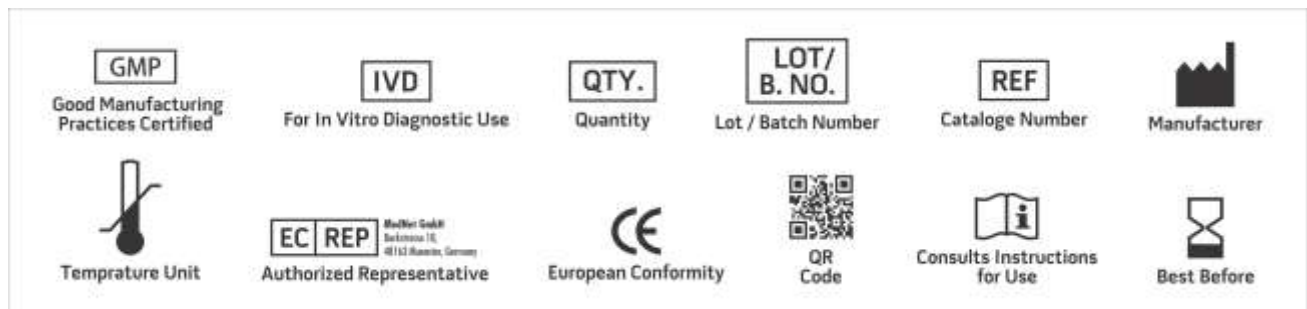

NOTE: Please consult the Material Safety Data Sheet for information regarding hazards and safe handling Practices.

***For Lab Use Only**
Revision: 28 Feb., 2022